

10-2 Details of the Function Items (For all models)

Function item	Parameter	Description				
<i>StYP</i>	<i>HtYPE</i> •	Portable type (This function can use: <i>H-Fnc</i> , <i>St-b</i> , <i>HLd-t</i> , <i>CP</i> , <i>bEP</i> , <i>CP H</i> , <i>CP Lo</i>) Factory setting of the SV-H series				
Change types	<i>RtYPE</i> •	Stand type (This function cannot use: <i>H-Fnc</i> , <i>St-b</i> , <i>HLd-t</i> , <i>CP</i> , <i>bEP</i> , <i>CP H</i> , <i>CP Lo</i>) Factory setting of the SV-A series				
<i>Cond</i>	0	Follows the viscosity changes quickly. (Prone to vibration)				
Condition	1 •	Follows the viscosity changes slowly. (Stable values)				
	2					
<i>Unit</i>	0 •	Viscosity	mPas	Temperature	°C	Factory setting: SV-1A/1H/10A/10H=0 SV-100A/100H=1
Unit upon power-on	1 •		Pas			
	2		cP			
	3		P			
	4		mPas			
	5		Pas			
	6		cP			
	7		P		°F	
<i>Pnt</i>	0 •	Dot		With "Comma" selected, the separator for CSV format will be ";" (semicolon).		
Decimal point	1	Comma				
<i>Fnc</i>	0 •	Switches viscosity units.				
[MODE] key function during measurement	1	Switches between the temperature display and the measurement elapsed time display.				
	2	Does not display the digits that are beyond the decimal point.				
<i>Prt</i>	0 •	Key mode		Press the [PRINT] key to output data.		
Data output mode	1	Auto print mode		Outputs automatically when the [STOP] key ends the measurement.		
	2	Stream mode		Continuous output during measurement. Outputs the viscosity only when D.P. format is selected.		
<i>TYPE</i>	0	A&D standard format		For AD-8121B MODES 1&2		
Data output format	1 •	D.P. format		For AD-8121B MODE 3		
	2	CSV format		For a personal computer		
	3	RsVisco format		For graphing program RsVisco		
<i>S-Rt</i>	0	No output		Available only for D.P. format		
Measurement elapsed time output	1 •	Output				
<i>S-td</i>	0	No output		Available only for D.P. and CSV formats		
Date/time output	1 •	Output				
<i>S-Ed</i>	0	No output				
Other output	1	Outputs remarks.		Available only for D.P. format		
	2 •	Outputs remarks, Device ID information and signature.				
	3	Outputs ID number.		Available only for CSV format		
<i>PUSE</i>	0	No pause				
Pause at data output	1 •	Pause (Approx. 2 seconds)				
<i>ErFnc</i>	0 •	Usually use this parameter.				
Reserved	1					
	7					
<i>H-Fnc</i>	0	Off				
Function of portable type	1 •	Auto hold function				
	2	Timer mode function				

• Factory setting